


Transferring Files from the Spectrometer PC

1. Open a file browser by double-clicking on the **Home** folder icon
2. Go to **File→Connect to Server...**
3. Enter the following into the window:

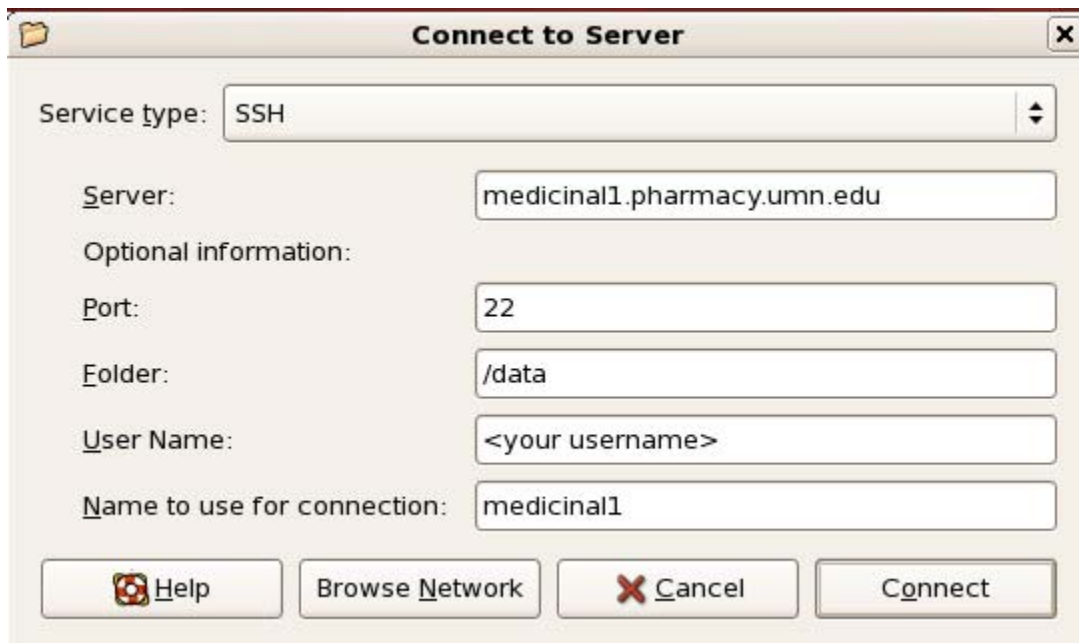

4. Click [**Connect**] and enter your password when prompted. A window should open showing the **/data** directory from the spectrometer computer. Files can be copied and pasted to the offline PC as needed.
5. After performing this procedure the first time, an icon will appear on your desktop. You can connect to the spectrometer PC in the future simply by double-clicking that icon and entering your password when prompted.

Note: The same procedure may be used to connect to ITDD. Change the following:

Server: itdd400.ahc.umn.edu

Folder: /opt/topspin/data

Name to use for connection: itdd400