

Activating BioPack in VNMRj 1.1D

1. Open the VNMRj classic interface of the software.
2. In the menu system, select [**Main Menu**] [**SetUp**] [**Activate Biopack**] [**Activate & copy vnmr probe file**].
3. This will start the BioPack removal macro that will back up any files from older versions of BioPack, Protein Pack or RNA pack. Another window will open and in that window at the ok to proceed prompt select [**y**]. The macro will check for any legacy files and back them up. After it is finished, hit the return key and the window will close.
4. At this point, the ghn_co parameter set should have loaded and it can be checked for accuracy. In the menu system, select [**Update Probe file with these Parameters**].
5. Wait for the activation process to complete.