

Use of Medicinal Chemistry PC Offline Computer

The intent of this computer is that it be used as a general access computer that supports one license for MestReC. MestReC is a third party processing software available through the Supercomputing Institute. The PC is connected via WinSCP to transfer files from the offline (medicinal2). The PC can also be used to back up files onto any flash drive and onto CD.

MestReC software

<http://www.mestrec.com/>

General Rules:

1. DO NOT LOGOUT OR SHUTDOWN THIS COMPUTER.
2. Keep all of your files in a single folder in “Documents” that is named with your name or username.
3. Please close all applications when you finish.
4. Do not store an excessive amount of data or files on this computer.
5. Files on this computer are not backed up by the facility. You are responsible for your own data.

Using WinSCP to transfer files:

1. Double click on the WinSCP icon.
2. Enter the hostname: medicinal2.pharmacy.umn.edu
3. Enter your login name and password for the offline Sun.
4. Login using the setting SFTP (allow SCP fallback).
5. This opens two windows – one for files on the PC and one for files on the Sun. Drag and drop the data you want to move from one window to the other.
6. Be sure to quit the connection when you are finished.