

Writing Data to a Floppy Disk on a Sun

The Ultra 10 (nmr1) in the Biochemistry NMR lab has a floppy drive that can be utilized for data storage of small files.

To use the floppy drive:

1. Insert the floppy disk into the drive.
2. Select the file manager icon from the CDE menu bar.
3. Choose "Check for Floppy".
4. A window for the floppy disk should open.
5. Then use the regular File Manager to drag and drop files to copy files to the floppy disk. The file manager can be accessed by clicking the right mouse button on the wallpaper and selecting "File Manager".
6. **Important!** When finished and want to eject the disk, go to the file window for the floppy and choose "File", "Eject". The message will about being able to manually eject the floppy now.
7. Use the eject button to remove the floppy disk.